

# At-a-glance Guide to the SenSci Data Logger

---

## Main Menu



**Meter Events**  
Displays data from  
Sensors but doesn't  
store results



**Log Events**  
Displays and  
stores data  
from sensors



**Time Events**  
Stopwatch feature  
and records times  
for digital sensors



**Data Files**  
View, analyse,  
upload and delete  
stored data files



**Settings**  
Adjust the  
system  
settings


---

## To Start Metering

1. Select



2. Select

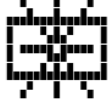

Snapshot

If you want to  
capture the data  
at that moment

or




Continuous

To get a continuous  
reading.

---

## To Start Logging

1. Select

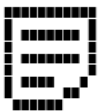

2. Select



Autolog

This logs at  
the rate set  
last time.

or



Manual Log

Logs whenever  
you hit the Select  
button

If you want to set the length of time you want to log for and how often you want the logger to take a reading.


1. Select



Parameters

to say how  
often the Logger  
should log

Select

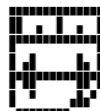

Lograte

This logs at  
the rate set  
last time.

Select




Next

to define the  
duration of the  
experiment.

---


## To upload data from the SenSci Data Logger

1. Select

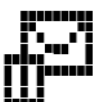


Data Files

2. Select

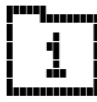

The  
Experiment  
you want to  
upload

3. Select

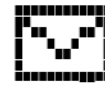

Upload